

deconex® 41 CLEAR

Acidic, ecological rinse aid

For use in washing machines



Application

deconex 41 CLEAR is used for rinsing laboratory glassware and labware in washers after the cleaning step. deconex 41 CLEAR is also used for reprocessing of medical instruments. However, this application is restricted to places where the quality of the water for the final rinse does not allow the use of a neutral rinse aid.

Properties

deconex 41 CLEAR neutralizes residues of alkaline solution and prevents lime deposits in the machine as well as water spots on washed items. Carefully selected surfactant compounds lend the product excellent wetting and clear rinsing properties.

In particular, glassware is provided with a nice shine. Moreover, the application of deconex 41 CLEAR accelerates the drying process of washed goods.

Ingredients

Organic acids, surfactants

Chemical/physical data

pH	1% solution in demineralized water	approx. 2.6
Density	concentrate	1.09 g/ml
Appearance	concentrate	transparent, light yellowish liquid
Titrate procedure	with 0.1 n potassium hydroxide	indicator: phenolphthalein
Consumption of 0.1 n potassium hydroxide	for 100 ml fresh washing fleet	35 ml/1% (w/w) of concentrate

Dosage

0.5-1.5 ml/l
depending on washing goods and on the quantity of alkaline residues

For reprocessing of medical instruments the dosage should be well controlled in order that the pH does not fall below 6.

Material compatibility

Suitable for:
Laboratory glassware, ceramics, stainless steel, aluminium, synthetic material

For materials not mentioned please make your own specific compatibility tests or consult Borer Chemie AG.

deconex® 41 CLEAR

Availability

Please ask your local representative about current container sizes.

Containers, screw caps and labels are made of recyclable polyethylene. Dosing aids such as pumps, dispenser, measuring jugs etc. are available. Please ask for corresponding documents.

Additional information

For information concerning safety at work, storage and waste disposal/effluent, please consult the corresponding safety data sheet.

Take advantage of our vast know-how! Please, contact us for further information regarding your specific application.



11924 Forest Hill Blvd., Suite 22-162, Wellington, FL 33414
Tel. +1 561 784 56 43, Fax. +1 561 880 68 97
www.racoma.com, contact@racoma.com

Manufacturer:

Borer Chemie AG

Gewerbestrasse 13, 4528 Zuchwil / Switzerland
Tel +41 32 686 56 00 Fax +41 32 686 56 90
office@borer.ch, www.borer.ch

All information provided is based on our current knowledge and it does not constitute a legally binding assurance of specific product properties.